

TechBriefs

Savannah River National Laboratory

U.S. DEPARTMENT OF ENERGY • SAVANNAH RIVER SITE • AIKEN • SC

srnl.doe.gov

At a glance

- > Collects microorganisms, chemical agents, radioactive particles
- > Collection efficiency of >90 percent
- > Collects particles of < 0.3 to 10 micron diameters
- > Maintains viability of live agents
- > 300 liters per minute flow rate
- > Consumes <12 watts of power
- > Uses a 12 volt battery
- > Measures 6" x 10" x 21"
- > U.S. Patent 6,955,075

Contact Information

Dale Haas

Commercialization Manager

Savannah River National Laboratory
Bldg. 773-41A, Room 238
Aiken, SC 29808

Phone: 803.725.4185

FAX: 803.725.4988

E-mail: dale.haas@srnl.doe.gov



Aerosol-to-Liquid Particle Extraction System (ALPES)

A new device uses electrostatic precipitation to collect and concentrate airborne agents in a liquid sample for onsite or laboratory analysis. The collection efficiency is 85 percent to 92 percent depending on the size of the particles. The Aerosol-to-Liquid Particle Extraction System (ALPES) is designed to collect chemical agents; radioactive particles; microorganisms such as spores, bacteria, and fungi; and molecules and other substances associated with explosives.

New device collects deadly agents for quick identification

Agents that could be used in chemical and biological warfare tend to disperse widely in the air when released. Quickly collecting a concentrated sample of these agents is critical to detecting them before they reach harmful dose levels in the air. Also, methods of detection such as by polymerase chain reaction are enhanced when the collected agents remain alive. Other collection devices using wet cyclone designs are larger and heavier, have higher power demands, and are noisy, thereby precluding unobtrusive uses. Devices using impactors for collection do not maintain the viability of live agents.

Uses proven particle separation technique

The new ALPES takes advantage of the long-proven use of wet electrostatic precipitation to separate particles from air. The device comprises an ionization section atop a tubular collection electrode enclosed in a column. A pump pulls air through the vertical column at a flow rate of up to 300 liters a minute.



TechBriefs

Savannah River National Laboratory

Technology transfer

The Savannah River National Laboratory (SRNL) is the U.S. Department of Energy's (DOE) applied research and development laboratory at the Savannah River Site (SRS).

With its wide spectrum and expertise in areas such as homeland security, hydrogen technology, materials, sensors, and environmental science, SRNL's cutting edge technology delivers high dividends to its customers.

The management and operating contractor for SRS and SRNL is Savannah River Nuclear Solutions, LLC. SRNS is responsible for transferring its technologies to the private sector so that these technologies may have the collateral benefit of enhancing U.S. economic competitiveness.

Recirculating liquid concentrates collected agents

A reservoir at the bottom of the column contains liquid that is pumped up through the inside of the collection tube, which is charged at 8,000 volts. The liquid flows over the top of the charged tube and down the sides, collecting the ionized particles from the outside surface of the tube. The flow rate maintains a continuous film of liquid on the outer surface of the tube. Constant recirculation of the liquid through the tube concentrates the collected particles.

A valved sample loop enables diversion of the liquid to an online analyzer or to a sample vial for transport to a laboratory for analysis.

Low power, quiet operation

With minimal pressure drop, the device consumes less than 12 watts of power and operates quietly.

Adaptable for selected analyses

The recirculating liquid can be customized for specific situations. A buffered saline solution or a nutrient solution will maintain biological agents in a viable state, enabling faster and more accurate analyses.

Partnering opportunities

SRNS invites interested companies with proven capabilities in this area of expertise to enter into a licensing agreement with SRNS to market this nuclear material detection system. Interested companies will be requested to submit a business plan setting forth company qualifications, strategies, activities, and milestones for commercializing this invention.

Qualifications should include past experience at bringing similar products to market, reasonable schedule for product launch, sufficient manufacturing capacity, established distribution networks, and evidence of sufficient financial resources for product development and launch.

SRNL-L9100-2008-00559



Savannah River National Laboratory[™]
OPERATED BY SAVANNAH RIVER NUCLEAR SOLUTIONS

The Savannah River Site and the Savannah River National Laboratory are owned by the U.S. Department of Energy, and are managed and operated by Savannah River Nuclear Solutions.